

Instructions for online submission of Shimadzu MALDI 8020/8030 spectra for comparison with the Mabriteccentral database

1. How to retrieve ascii txt files

All spectra are stored in the export folder (Contact Shimadzu fur further details).

2. How to upload ascii txt files from shimadzu to mabriteccentral.com

😵 Sign In with Auth0 🛛 🗙 🖸	Google x +	- C ×
(←) → C ^a (a)	🛈 🔒 https://www.google.com//webhp?hl=en&sa=X&vved=0ahUKEwj4w53ojcXiAhUK-qQKHeCvAcgQPAgH 🛛 😶 😨 🟠	In © ≋ ≡
	Coogle Search I m Feeling Lucky Google offered in: Deutsch Français Italiano Rumartach	Grnail Images 🔀 💙
	Open Login via <u>www.mabriteccentral.com</u>]
Switzerland		
Advertising Business About		Privacy Terms Settings

2.1 Access mabriteccentral.com via browser

2.2 Enter your credentials



$\leftarrow \ \ \rightarrow \ \ G$	○ △ ē ² https://app.mabriteccentral.com//ogin	☆	© ± 0% £
	mabritec CENTRAL		
	Welcome log in to continue to mabriteccentral.com		
	Ernail address Password Forgot password?		
	Continue		
	Enter username and password		
	Price and License Legal notice Privacy Terms of Service mabritec - gesellschaft für massenbasierte rasche identifikation	nstechnologien	

2.3 Upload your mzml or ascii files

CENTRAL	① Identification ① Mosquito Vectors ① Spectra Quality 🗮 Orders 🛄 Tutorial 옫 Account 더 Logout
	Identification Identify your unknown spectra or confirm your Identification result. Shimadzu 8020 / 8030 (ascii)
	01. Select your machine type: O Bruker Biotyper O Shimadzu 8020 / 8030 (ascir) O Clover BioSoft O Vitek MS (mzml) O Vitek MS Prime (ascii)
	02. Select a database: 03. Add customer referen Choose database all bacteria
	O4. Upload files:
	Drag and Drop files here, or click to select them from your computer or select your bruker compass experiment <u>her</u> (file size must not exceed 30MB)
	Clear Send Move the ascii files to be transmitted into the "Drag and Drop" field. Or navigate to the corresponding folder by "clicking" on the "Drag and Drop" field and select all ascii files.

mabritec CENTRAL

2.4 Select/Deselect

$\leftarrow \ \rightarrow \ G$	O 🔒 🔤 https://app.mabriteccentra	.com/upload	다 ⓒ 노 💿 한
		01. Select your machine type: O Bruker Biotyper O Vitek MS (mzml) O Vitek MS (ascii)	
		02. Select a database:	
		All Bacteria ~	
		03. Upload File:	tions
		Drag and Drop files here, or click to select them from your computer (file size must not exceed 10MB)	2
	1	REFERENZ-D11.txt 4.87 kB ×	
		REFERENZ-D12.txt 4.86 kB ×	
		Clear Send 3	
	1. The selec	ted files are displayed.	
	2. Files that	were selected by mistake can be dele	eted by means of "delete".
	3. Otherwis	e launch upload and identification by	pressing "send".

2.5 Results / PDF and CSV export

→ C nabritec ENTRAL	A # http://app.mabriteccentral.com/order/5615	4c7e-b041-4at7-892a-f7f9d5a9a651		ti ① Upload 듣 Orders ∷	⊖ ⊻ w
	Order ID: 56114c7e-b041-4af7-892a-f7f9d5a9a651 Order date: 20.3.2023, 11:33:24 Number of spectras uploaded: 8 Database used: All		CSV export		
	Filename	Sample Name	Qualified result	Non qualified result	
	042_Mabritec_AXIMA_1.09.mabr_E QA20_00022.7.txt more details	042_Mabritec_AXIMA_1.09.mabr_E QA20_0002.2J1	Escherichia coli	0	
	042_Mabritec_AXIMA_L09.mabr_E GA20_00022J2.txt more details	042_Mabritec_AXIMA_1.09.mabr_E QA20_0002.2J2	• Escherichia coli		
	Results are displaye	ed as a list and ca	in be exporte	d as a PDF or a CSV f	file

Mabriteccentral - mabritec AG - Lörracherstrasse-50 - 4125-Riehen - Switzerland www.mabriteccentral.com